

“Blue” Physical Property Measurement System (PPMS) – General Instructions & Info

All users must receive training for the PPMS and for whichever specific options they wish to use. Reserve all necessary time in FOM and log on and off when starting and finishing your experiment.

This document contains general system information and control instructions. The PPMS Hardware manual and the individual option manuals contain more detailed information. Please refer to the specific option SOPs for information on changing options/samples and making measurements. Please consult the manuals and/or the facility staff for help and troubleshooting information.

Model 6000 Controller

This control unit sits in the PPMS cabinet and is the main controller for general hardware operations. While system controls are generally performed via the MultiVu software, all hardware control operations can also be performed using the Model 6000. This is primarily useful if the computer/software crashes or is unavailable for any reason. In some cases, changing options requires interchanging cable connectors on the back of the Model 6000. See the specific option SOPs for details on such changes.

GasMon32

While not typically used for normal system operations, this program can be useful for troubleshooting issues with the gas/cryo system or for determining the current state of a specific system element. It also serves as a good schematic of the chamber and helium dewar. To open, click the Gasmon32 taskbar icon (yellow “32”). In addition to the status of various valves, the program includes read-outs for the helium level and all system pressure and temperature sensors.

MultiVu

Main hardware controls

PPMS MultiVu is the main interface system for all hardware and measurement controls. The bottom of the window (see below) contains the main hardware system indicators, which can be used to access their respective controls by clicking the appropriate area. From right to left:

Sequence Run	35.994 K, Near	100.59 Oe, Persistent	12.24 Torr
Seq: ACDC-Tsweep.seq	Set: 36.000 K	Set: 100.0 Oe	Purged
3: Scan Temperature	2.00 K/min, Fast Settle	100.00 Oe/sec, Linear	63.10% He

- **Chamber** – Indicates the sample chamber pressure, the gas control state, and the helium level. Normal operating pressure is 7-10 torr (helium). Control buttons include *Seal* (seal off chamber from pump and gas inlets), *Purge/Seal* (chamber is purged with helium gas and pumped three times, then sealed), *Vent/Seal* (chamber is purged with helium gas, then sealed), *Pump Cont.* (chamber is pumped continuously), *Vent Cont.* (chamber is continuously flooded with helium), and *HiVac* (chamber is pumped with cryopump – see below).
- **Magnet** – Indicates the magnetic field and state of the magnet, the field setpoint, field ramp rate and approach mode. Control fields include: Setpoint (-140,000 to 140,000 in Oe where 1 T = 10,000 Oe), Rate (0 to ~130 Oe/sec), Approach (Linear/No O’shoot/Oscillate – see manual for definitions), and Mode (Persistent/Driven). The normal operating state is Persistent, meaning a current (or no current for 0 field) is flowing through the superconducting electromagnet without any driving voltage. When the field is changed, a switch is activated

(the state briefly reads “SW Warm”), a voltage is applied to change the drive current and thus the field (the changing field is indicated along with “Ramping”), then when the field is reached, the switch is opened (“SW Cool”) and the field stabilizes. Under Driven mode, the switch always remains closed. This allows for faster changing between fields, but uses significantly more helium due to ohmic heating of the active switch. Regardless of the switch mode used, there is a brief delay while communication is established with the external power supply, and the state is “Undefined”. Persistent mode should be used to conserve helium; please consult with facility staff if you feel that Driven mode is necessary for your measurement.

- **Temperature** – Indicates the system temperature and control state, the temperature setpoint, and the ramp rate and approach mode. Control fields include Setpoint (1.9 to 400 K), Rate (0 to 20 K/min), and Mode (Fast Settle/No O’shoot – see manual for definitions). Note that the software will accept ramp rates up to 20 K/min to be input, but the system will likely not actually ramp this fast. Actual maximum ramp rates depend on the option being used and the temperature range. The temperature should not be immediately changed from high temperatures to < 10 K. In such cases, set the temperature to 10-15 K, wait for 10-30 minutes, then continue to the lower temperature. This practice should also be used within sequences.
- **Sequence** – See below.

Data and sequences

While hardware operations can be controlled directly through the Model 6000, MultiVu is required for saving data and running automated sequences. In general, data from any measurement option can be saved for immediate measurements (these are generally handled through the option control windows – see option SOPs for details) or for measurements as part of a sequence. In general, the user has to specify to have data from immediate measurements saved, and data from measurements in sequences is automatically saved. Data is saved in .dat files, which are comma-delimited text files that can be imported into any standard analysis program. Opening a .dat file within MultiVu (this can be done in the main software window, or by clicking the **View** button within the data file area of the active option control window) will open a graph by default. Right clicking brings up a menu which includes Data Selection for changing the data axes and ranges (multiple Y axes are available) and Table View for viewing the data in table form. Once a file has been opened as a graph, MultiVu also creates an additional file with .gph appended to the file name; this file automatically saves the graph with the most recent data and axes.

Sequences are sets of control and measurement commands which can be used to run automated experiments. They are saved as .seq files, which appear in MultiVu as a simple series of command lines. The right side of the MultiVu screen is the menu of sequence commands. Double-clicking any item will add that command to the sequence above the line in the sequence which is highlighted. Double-clicking on a line within a sequence allows the command parameters to be edited. All the operations listed at the top of this page and many more are available. Commonly-used commands include temperature and field scans, during which measurements can be taken at specified intervals. The list of available measurement commands depends on the active option, and generally needs to be expanded by clicking “+” next to the option name. After writing a sequence, it must be saved, at which point clicking **Run** on the left side of the screen or the “play” button on the top toolbar will start the sequence. The current sequence status is given in the bottom (see above) and left sections of the screen. Some notes on sequences:

- Individual or multiple (using Ctrl-click or Shift-click) lines can be selected and cut, copied, pasted, or deleted.
- Right-clicking a selected line or set of lines allows it to be disabled, so it can be excluded from a running of the sequence without being permanently deleted from it.
- All temperature and field scans will automatically have an “End scan” line at the end. Any lines within the scan will be indented and will be executed at each point within the scan.

- Within a temperature or field scan, Sweep is an additional Mode along with the two or three outlined above. In this mode, measurements are made as the specified temperature or field is reached, and the temperature or field continues sweeping without ever settling; ramp rates should be selected with the duration of the individual measurements in mind in order to avoid a significant T or H change within one measurement.
- At the end of all sequences, set the system temperature to 300 K and the magnetic field to zero Oe (persistent).

Cryogen System

The PPMS is cooled by a liquid helium dewar with a liquid nitrogen jacket. Boil-off from the helium is collected by the reliquefier and returned to the dewar. The nitrogen vents into the room through pressure relief valves. The valve on the left side of the dewar lets off exhaust at 1 psi and is frequently active. Users should use the hot air gun to remove excess ice, and in particular, make sure that ice does not freeze the valve open (air can enter into the dewar) or closed (dewar can over-pressurize). The helium level is maintained via periodic fills run by the facility staff and should always be above 70%, and ideally above 80%; please notify the staff immediately if you observe a lower level.

Cryopump

The cryopump can be used to reduce the sample chamber pressure to $\sim 10^{-5}$ torr. This can be used for any measurement where a vacuum environment is preferable (keep in mind that the system normally relies on the helium gas in the chamber for thermal exchange) and is required for use of the Thermal Transport Option (TTO) or Helium-3 option. The cryopump manual includes details on the theory of operation and the physical configuration of the pump, including detailed operating instructions in chapter 3. If the chamber has been opened, a Purge & Seal should be performed as usual, and then HiVac can be initiated using the chamber controls. The system will initially enter “Pre-HiVac” mode for 1-2 minutes, then switch to HiVac mode, after which the base pressure should be reached within a few minutes. For puck-based measurements at low temperature using the cryopump, the contact baffle (stored in cryopump user kit) should be attached to the end of the baffle rod. If the chamber fails to reach HiVac status within ~ 10 minutes, please notify the staff; the cryo pump may need to be regenerated.

Physical Property Measurement System (PPMS) – Resistivity (DC)

All users must receive specific training for the Resistivity option regardless of training on other options. Reserve all necessary time in FOM and log on and off when starting and finishing your experiment.

The Resistivity option is used for making basic DC resistance measurements by applying a current and measuring the resulting voltage. Resistance measurements can also be made using the AC Transport option. Which option is more appropriate depends on several factors, including whether an AC signal is desirable/acceptable, the resistance of the device(s) to be tested, and the number of devices to be tested. Consult the system documentation and/or facility staff to help determine which option is appropriate. The three “Transport...” PDF documents on the Red PPMS computer desktop are a particularly useful resource for making electronic transport measurements.

Sample Preparation

The Resistivity option performs a 4-point resistance measurement on up to 3 channels at a time. The standard Resistivity puck (see right) includes pre-mapped pads for + and – current and voltage for Channels 1 – 3. The recessed area measures 11.6 mm x 13.6 mm (with rounded corners), so samples that fit within this area will work best. The base is conducting, so an insulating layer such as Kapton tape will be necessary for samples with conductive back sides.

Electrical contact between the sample and the desired channel input pads can be made via wire bonding, soldered or silver-painted wires, spring-loaded press contacts (consult staff on use of these), or other methods. Alternatively, samples can be mounted on the universal puck, which has a blank face and no pre-configured channel set-up. This can be useful for larger samples or if no Resistivity pucks are available. The pin-to-channel



correspondence for the user puck is posted on the bulletin board. The Resistivity or universal puck can be used in combination with the P150 User Bridge (the red face plate corresponds to the Resistivity option channels) to test sample-to-puck connections in advance of insertion into the PPMS using a hand-held ohm meter or other device.

Activation

1. In the MultiVu software, de-activate any other active option, and physically remove the cables/hardware as necessary according to the respective SOP.
2. Attach the Resistivity LEMO connector to the back of the probe head.
3. Attach the other end of the cable to the P1 User Bridge port on the Model 6000.
4. Click **Utilities – Activate Option...** and select and activate the Resistivity option. The Resistivity Option control window should appear if the option has been successfully activated.
5. If not immediately inserting a sample, insert the baffle/flange rod and clamp it in place, then purge and seal the system (if necessary). Or, go to step 3 under Measurement Procedure to install a sample.

Measurement Procedure

Refer to Chapter 3 of the Resistivity manual for information on performing measurements. Once the option is activated, the Resistivity Option control window contains most commands needed to set up and perform a measurement. To insert a sample:

1. Click the **Install/Remove...** button and follow the on-screen instructions.
2. When the system is vented, remove the baffle/flange rod.
3. Visually verify that there is not a puck in the system, and if there is, remove it.
4. Place the puck in the puck insertion tool:
 - a. Flip the switch on the handle up to open the ring.

- b. Put the puck into the bottom of the tool, sample side up.
 - c. Hold the puck in place while flipping the switch down to lock the puck in place.
5. Using the tool, slowly lower the puck into the probe. When it reaches the bottom, rotate it until the key aligns with the key slot (this happens with the key facing forward), then press down to engage the pins. Do not press too hard – if the pins are not engaging properly, there may be damage which could be worsened.
6. Flip the switch on the tool up and withdraw the tool.
7. Replace the baffle/flange rod and finish the install procedure in the software to purge and seal the system.
8. Place the puck insertion tool back in the tube in the storage unit.
9. Under the Datafile header, click `Browse...` and set up a file location and name. Once this is defined, you will be prompted to enter sample info for any channels to be used. This includes the length and cross-sectional area to be used in the resistivity calculation; these can be ignored if the sample geometry is not appropriate for a simple bar geometry calculation of resistivity.

When the sample is inserted, real-time resistance readings can be seen via the bridge set-up, which must be defined by clicking the `Bridge Setup...` button. In the subsequent window, parameters for each of the three channels (plus Channel 4, which is not used) can be specified. See page 2-10 of the Resistivity manual for details. Note that Channel Set dictates whether the excitation current is being applied or not, and that the measurement can be limited by any one of current, power, or voltage depending on the specific limits specified and the sample characteristics (the graph on page 2-12 provides a nice illustration of this). Despite the “DC” designation, all measurement excitations are actually single-pulse measurements at 60 Hz (corresponding to the input line frequency), and Drive Mode dictates whether this is a single positive pulse (DC) or a full-cycle bi-polar pulse (AC, which is the default). The right side of the window gives the “Status” of each channel – the specified parameters are applied whenever the channel is on, and the current and resistance are displayed.

For data to be saved, a “measurement” (as opposed to a simple real-time indication) must be made, either in immediate mode or as part of a sequence.

- Immediate measurements can be made by clicking the `Measure...` button and using the Resistivity Quick Measure window. The channel(s) to be measured and the number of readings to average are specified. Note that the bridge can be re-defined within the measurement window by clicking `Bridge Setup...` Click `Measure` to begin the measurement. Results are displayed and can be saved to the specified file (see above) by clicking `Save`.
- For information on setting up and running sequences, refer to the PPMS general instructions. Resistivity measurement commands are listed under “Resistivity” (click the + to expand the list of choices) and are similar to immediate measurements. The Resistivity measurement definition window again allows the bridge set-up to be defined using the Bridge Configuration tab, and the Data Fields tab allows the user to select what parameters and results (from both the measurement and the system hardware) will be saved to the file.

To change samples, set the system temperature to 300 K and the magnetic field to 0 Oe (persistent); this should be done at the end of each sequence, but can be done manually if necessary. Use the sample change sequence as described above (with slight differences for puck removal). When finished, the Resistivity option can be left active in the system.

De-activation and Removal

1. Click `Utilities - Activate Option...` and select and de-activate the Resistivity option.
2. Remove the Resistivity LEMO connector from the back of the probe and hang the LEMO connector from one of the hooks on the wall.


Physical Property Measurement System (PPMS) – AC Transport (ACT)

All users must receive specific training for the AC Transport option regardless of training on other options. Reserve all necessary time in FOM and log on and off when starting and finishing your experiment.

The AC transport option can be used for making several different measurements with AC excitation: resistivity, I-V curves, Hall Effect, and critical current. Resistance and (implicitly) Hall measurements can also be made using the Resistivity option. Which option is more appropriate depends on several factors, including whether an AC signal is desirable/acceptable, the resistance of the device(s) to be tested, and the number of devices to be tested. Consult the system documentation and/or facility staff to help determine which option is appropriate. The three “transport...” PDF documents on the Red PPMS computer desktop are a particularly useful resource for making electronic transport measurements.

Sample Preparation

The AC Transport option performs measurements using 4 or 5 contact points on up to 2 channels at a time. The standard ACT puck (see right) includes pre-mapped pads I+, I-, V_A+, V_B+, and V- on both Channels 1 & 2. The recessed area measures 10.8 mm square (with rounded corners), so samples that fit within this area will work best. The base is conducting, so an insulating layer such as Kapton tape will be necessary for samples with conductive back sides. Electrical contact between the sample and the desired channel input pads can be made via wire bonding, soldered or silver-painted wires, spring-loaded press contacts (consult staff on use of these), or other methods. Alternatively, samples can be mounted on the user puck, which has a blank face and no pre-configured channel set-up. This can be useful for larger samples or if no ACT pucks are available. The pin-to-channel correspondence for the user puck is posted on the bulletin board. The ACT or user puck can be used in combination with the User Bridge (the blue face plate corresponds to the ACT option channels) to test sample-to-puck connections in advance of insertion into the PPMS using a hand-held ohm meter or other device.

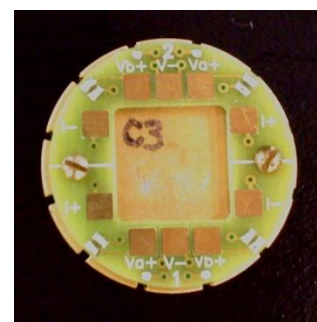

Activation

1. In the MultiVu software, de-activate any other active option, and physically remove the cables/hardware as necessary according to the respective SOP.
2. Click *Utilities – Activate Option...* and open the ACT connection diagram and connect all cables accordingly. Note that if the ACMS option has been used more recently than the AC Transport option (check the log book to determine this if necessary), two cables on the back of the Model 6000 controller will need to be changed.
3. Select and activate the AC Transport option. ActStatus.log will pop up to indicate the activation status.
4. If not immediately inserting a sample, insert the baffle/flange rod and clamp it in place, then purge and seal the system (note – this is not necessary if switching from the Resistivity option). Or, go to step 3 under Measurement Procedure to install a sample.

Measurement Procedure

Refer to Chapter 5 of the AC Transport manual for information on performing measurements. Once the option is activated, the tabbed AC Transport Control Center contains most commands needed to set up and perform a measurement. To insert a sample:

1. On the Install tab, click the *Vent* button and follow the on-screen instructions.
2. When the system is vented, remove the baffle/flange rod.

3. Visually verify that there is not a puck in the system, and if there is, remove it.
4. To place the puck in the puck insertion tool:
 - a. Flip the switch on the handle up to open the ring.
 - b. Put the puck into the bottom of the tool, sample side up.
 - c. Hold the puck in place while flipping the switch down to lock the puck in place.
5. Using the tool, slowly lower the puck into the probe. When it reaches the bottom, rotate it until the key aligns with the key slot (this happens with the key facing forward), then press down to engage the pins.
6. Flip the switch on the tool up and withdraw the tool.
7. Replace the baffle/flange rod and finish the install procedure in the software to purge and seal the system.
8. Place the puck insertion tool back on the wire shelving unit.
9. On the Data File tab, click `Browse` and set up a file location and name. Then, you will be prompted to enter sample info on the Sample 1 and Sample 2 tabs; click `Next >>` and `Finish` after each one, respectively.

All measurements made using AC Transport fall under one of four modes corresponding to the four buttons which are always visible at the bottom of the AC Transport Control Center (regardless of which tab is open): `Hall Coefficient`, `Resistivity`, `IV Curve`, or `Critical Current`. Clicking one of these buttons will open a new window which serves as the interface for that measurement. Since each interface and measurement procedure is unique, users should refer to the relevant sub-section of section 5.3 of the AC Transport manual for specific information on a measurement type. In general:

- Immediate measurements can be made by clicking the `Measure` button at the bottom of the measurement window. Results are displayed and can be saved to the specified file (see above) by clicking `Save`.
- Within the AC Transport Control Center, the Waveforms tab displays the expected and actual data waveforms for every measurement. This can be useful for seeing noise levels and determining if the input parameters and measurement range are appropriate.
- For information on setting up and running sequences, refer to the PPMS general instructions. ACT measurement commands are listed under “AC Transport” (click the + to expand the list of choices). When a measurement command is selected, the measurement interface includes tabs for Sample 1 and Sample 2; the measurement can be made on either or both samples (each with their own separate parameters) by clicking the respective Measure check-boxes, whereas one or the other is selected in an immediate measurement. The PPMS Data tab allows the user to select what parameters and results (from both the measurement and the system hardware) will be saved to the file.

To change samples, set the system temperature to 300 K and the magnetic field to 0 Oe (persistent); this should be done at the end of each sequence, but can be done manually if necessary. Use the sample change sequence as described above (with slight differences for puck removal). When finished, the AC Transport option can be left active in the system.

De-activation and Removal


1. Click `Utilities - Activate Option...` and select and de-activate the AC Transport option.
2. Remove the AC Transport LEMO connector from the back of the probe and, if necessary, the connectors from the back of the Model 6000 controller. Hang the LEMO connector from one of the hooks on the wall.

Physical Property Measurement System (PPMS) – Horizontal Rotator (HR)

The Horizontal Rotator is not technically an option; it is an accessory which can be used in conjunction with the Resistivity and AC Transport options. Any measurement that can be made with those options can be made with the rotator, with two exceptions: the current limit is reduced from 2 A to 0.5 A in ACT, and only two channels are available in Resistivity. The Horizontal Rotator consists of a motor and the rotator probe body. The motor control is always active, so it simply needs to be connected via the control cable, mounted on the system, and coupled to the rotator body (which is installed in the probe space) in order to be used. The probe can also be rotated manually outside the system using the motor coupling. The positions are defined such that the sample faces up at 0° and forward at 90°. Slightly more than one full rotation is possible so that the end positions are defined as -10° (going past 0 to 350 on the dial) and 370° (going past 0 to 10). An on-board thermometer is used rather than the main PPMS system thermometers.

Sample Preparation

The Horizontal Rotator uses special sample holders rather than the pucks which are used with the Resistivity or AC Transport options. Samples can be connected to them in the same manner as the pucks, and connections can be tested in the user bridge via the appropriate adapter (all parts referred to in this section are kept in the Horizontal Rotator accessory kit). The sample holders are shown below.

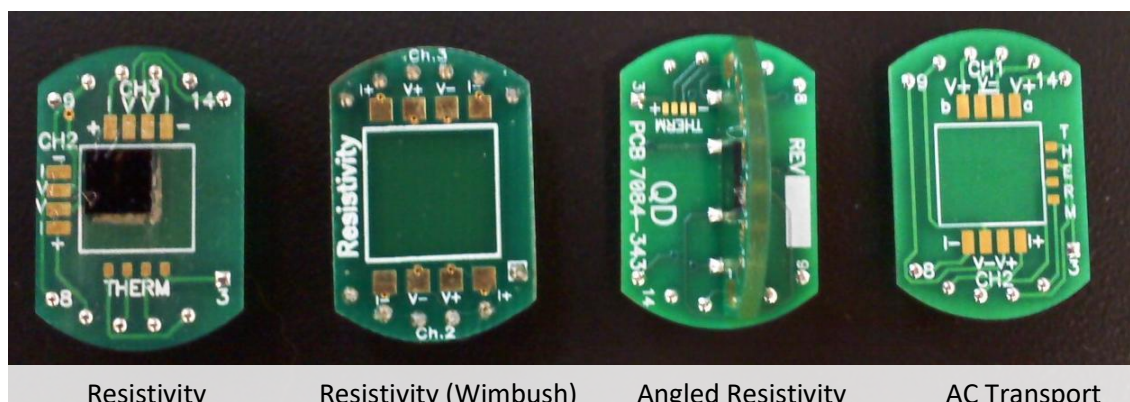


Resistivity – Channel 1 is designated for the thermometer – DO NOT USE THESE PADS. Channels 2 and 3 are for 4-point resistance measurements.

Resistivity (Wimbush) – Third-party version of the Resistivity holder with several improvements: larger sample area, clear “Resistivity” label, larger contact pads, and no exposed thermometer pads.

Angled Resistivity – Identical connections to the Resistivity holder, but with a second perpendicular board coming out of the main board so that the sample can be rotated within its own plane.

AC Transport – For 5-point Hall measurements, Channel 1 must be used; for other measurements, either Channel 1 or 2 may be used. Only the voltage pads are associated with the specific channels, and there are only two current leads which are shared by the channels and used for the active measurement at any given time. See page 7-4 of the AC Transport manual for details.

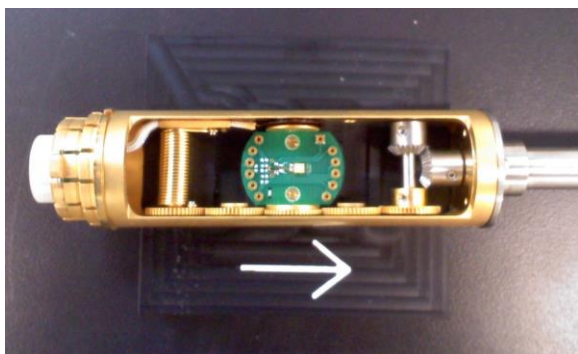

Once the sample is mounted to the holder, the holder must be attached to the rotator. Manually set the rotator position to 90° and use the rotator support tool to securely rest it on a flat surface as shown to the left. Then, press the sample holder into the rotator to engage the pins; note that the pin configuration is asymmetric, so it will only attach in one orientation. The sample removal tool is designed for this purpose and can grip the sample holders at the curved edges.

Rotator Insertion and Motor Set-up

1. Click `Utilities - Activate Option...` and open the connection diagram for the appropriate rotator/option combination and connect all cables accordingly.
2. If necessary, select and activate the Resistivity or AC Transport option.
3. If necessary, vent the system and remove the flange/baffle rod assembly.
4. Visually verify that there is not a puck in the system, and if there is, remove it.
5. Place the rotator motor assembly on top of the system abut do not secure the flange. Move the motor off to one side as shown to the right.
6. Turn the rotator to -10° .
7. Carefully insert the rotator into the probe space. When it is nearly in, lift the rotator support so that the rotator passes through it. Then, turn the full rotator body to the proper position (key/white line forward) and firmly push it down to engage the pins.
8. Push the support down to form a seal with the o-ring, then secure the flange in place by pushing the wings in.
9. Next, the rotator's on-board thermometer must be activated. Open the program PPMS Rom Config 32 (taskbar icon with magnifying glass and chip) and click the "Diag. all" check box. Then, click `Send to PPMS - Send Config` and open the appropriate configuration file (HR-HiRes-442 for Resistivity or HRACT-HiRes-442 for ACT). The User temperature on the Model 6000 will now be active, and the temperature in MultiVu will match it.
10. Open the motor control window by clicking `Instrument - Motion...`. The full range of -10° to 370° should be indicated on the scale, as well as the current motor position. Click `Go to Index` to move the motor to the end position of -10° .
11. The motor and the rotator should now be in the same position and can be coupled by moving the motor to the center and lowering it onto the coupler.



Rotator Control

The control window described in step 8 above is used for immediate control of the rotator. To move to an angle, simply type the desired angle in the "Move to:" box and click `Set`. At this point, an immediate measurement can be made using the active option as described in the appropriate option procedures. Note that there is a slight hysteresis, so a given position may differ by a few degrees depending on from which direction it is approached. **Do not use "Redefine Current Position:"**. If you think something is wrong with the position definitions, consult the facility staff. When using sequences, "Set Position" and "Scan Position" are available in the command menu; these are used similarly to the equivalent commands for temperature and magnetic field.

De-activation and Removal

1. First, the on-board thermometer must be de-activated. Open the program Mon6000 (taskbar icon with a blue 32) and enter the command line "usertemp 0". The User temperature on the Model 6000 will now be blank.
2. Vent and seal the system.
3. Lift the motor to de-couple it from the rotator and move it to the side.
4. Carefully remove the rotator from the probe space, turn it to 90° , and set it on the rotator support tool on the table.
5. Remove the motor assembly from the flange and place it securely on the storage rack.
6. Insert the flange/baffle rod assembly and purge and seal the system.
7. If necessary, de-activate and disconnect the active measurement option as described in the option procedures.
8. Remove your sample from the rotator and leave the rotator in the storage tube on the rack.


Physical Property Measurement System (PPMS) – Helium-3 Refrigerator (He-3)

All users must receive specific training for the He-3 option regardless of training on other options. Reserve all necessary time in FOM and log on and off when starting and finishing your experiment. All users must fill out the He-3 usage log in order to track the remaining He-3 gas pressure over time and to determine whether cryo-cleaning of the He-3 is necessary before starting.

The He-3 option allows sample cooling to $\lesssim 400$ mK. The sample is mounted on a dedicated circuit board which attaches to the He-3 insert in order to provide direct thermal contact with a closed-loop He-3 circuit. At low temperatures (below 20 K), the He-3 option is operated under high vacuum (via the cryo pump) in order to eliminate convective heat transfer; hi-vac mode is initiated automatically at the appropriate time. For details on the temperature control methods (which vary with temperature range) and operations, please consult the Quantum Design manuals. Note that our system uses the GPIB version of the option (as opposed to the CAN version). The upper temperature limit for this option is 350 K.

Sample Preparation

The drawing to the right shows the He-3 sample holder for transport measurements. Samples up to 1 square cm can be accommodated. All samples should be attached using low-temperature varnish; Apiezon grease may suffice for light samples. The sets of four contact pads correspond to Channel 1 (pins 3-6) and Channel 2 (pins 7-10) of the Resistivity measurement option.

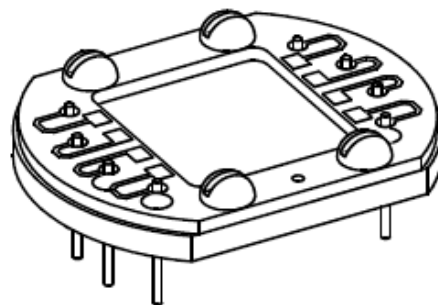

Set-up and Activation

1. Roll the cart up to the dewar and remove the cover. It needs to be close for the cable to reach. Remove the protective tube from the probe and the red cover from the o-ring.
2. Mount the sample holder onto the probe. Use Apiezon N grease to enhance thermal contact; if you plan to exceed 320 K, use Apiezon H. Using two short 0-80 screws and a 0.050" Allen wrench or ball driver, secure the holder in place; the drawing above shows four brass screws, but there will only be two, with two holes available for this purpose. [Screws, an Allen wrench, and small tubes of both greases are stored in the He-3 kit.]
3. If necessary, plug the power cord into the He-3 controller. It uses 220 V power, so be sure to use the dedicated power cord coming from the PPMS power distribution unit (attached to left side of tower) rather than a standard cord plugged into a wall outlet.
4. Make sure the green TANK supply valve is open.
5. If necessary, cryo-clean the helium-3 using the liquid nitrogen trap. Cleaning is required every two weeks if the option is regularly in use (facility staff will perform these when possible), or if the option has been idle for more than two weeks.
 - a. Fill the liquid nitrogen bucket using the low-pressure supply dewar in H21.
 - b. Slowly lower the cold trap into the liquid nitrogen.
 - c. Run the He3GasMon utility (located in C – QdPpms – He3Option) and launch the “Clean He3 with LN2 Trap” wizard and follow the on-screen prompts.
 - d. When directed to connect the trap, remove the protective caps from both fittings (trap and module) and apply a thin coating of Apiezon M grease to each face. Attach the fittings and then open the green CRYOCLEAN valve (note that the wizard will not prompt you to open the valve).
 - e. Proceed with the wizard and allow the cleaning process to complete (~20 minutes). When finished, close the CRYOCLEAN valve (again, the wizard will not say this) before disconnecting the trap.
 - f. Close He3GasMon.
6. If necessary, apply a thin coating of Apiezon N or H grease to the contact fingers along the probe body.

7. Vent and seal the PPMS and remove the flange/baffle rod.
8. Remove the probe from the cart and carefully lower it into the PPMS chamber with the label facing forward. Be careful to avoid putting too much tension or torque on the cable and watch out for the pipes above the chamber. For most users, it will be necessary to stand on the step stool. The probe should engage on the pins if aligned properly; otherwise, the brackets will not fit around the top of the chamber. You should feel the springs compress as the probe bottoms out. See photo for proper installed position.
9. Purge and seal the chamber.
10. Activate the He-3 option using the Option Manager. This will launch a control module which will list the initialization stages; once initialization completes, it will say that the option is controlling temperature. Note that the He-3 thermometer will occupy the User line on the Model 6000 and MultiVu (similar to other options with dedicated thermometers) and it will also occupy the Temp line on the Model 6000 (normally maintained by the system thermometers).



Measurements

While AC Transport measurements are possible using the He-3 option, at this point, we are only equipped for Resistivity measurements. Once the He-3 option is configured and activated per the above section, the Resistivity measurement option can be operated as it normally would be with a sample on a puck. Consult the Resistivity option SOP for details. Please note that the He-3 thermometer uses Channel 3, so if you want to run an existing sequence, all measurement steps need to have Channel 3 set to No Action rather than Off.

De-activation and Removal

1. De-activate the He-3 option using the Option Manager.
2. If the He-3 option will not be used in the next two weeks, the gas should be returned to the tank for storage using the “Clean and Secure He3 Gas” wizard in He3GasMon. Follow the instructions outlined by the wizard and record the He-3 pressure in the tank after completion.
3. Warm the system to 310 K and leave it there for at least 10 minutes.
4. Vent and seal the system.
5. Carefully remove the probe from the chamber and place it securely on the cart.
6. Cover the o-ring with the red cover and the probe with the protective tube.
7. Place the cover back on the cart and roll it back into the corner.
8. Put the flange/baffle rod back in the chamber, purge and seal, and return the system to 300 K.